

10. Setup



192.168.173.1:8085/admin

WiFi Access Point

Camera
Live image size: 640x480

Admin
 Require PIN
Set a PIN to restrict access to this page.

Networking
 Built-in WiFi Access Point
Password: 00000000
 Join existing WiFi

www.leica-microsystems.com



Local WiFi

Require PIN
Set a PIN to restrict access to this page.

Networking
 Built-in WiFi Access Point
 Join existing WiFi
Name (SSID):
Password:

AP: The HUB can act as a WiFi Access Point allowing the iPad connect directly - no infrastructure required. Though you will need to connect your iPad to the HUBs own WiFi network (DMshare-PT0001) which will prevent internet access.

Existing Network: you can connect the DMshare Hub to your existing WiFi network - the one your iPad is already connected to - so you don't need to switch back and forth for internet access.

Living up to Life




Leica DMshare Quick Start Guide


ryf ag  Ryf AG
Bettlachstrasse 2
2540 Grenchen
tel 032 654 21 00
fax 032 654 21 09
www.ryfag.ch

Order no.: English 934 020 Subject to modifications.
Copyright © by Leica Microsystems CMS GmbH, Wetzlar, Germany, 05/2014.
LEICA and the Leica Logo are registered trademarks of Leica Microsystems IR GmbH.


1. Data Transfer Hub

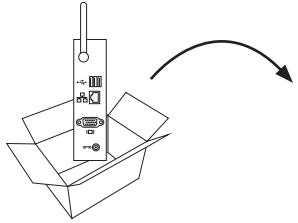

2. USB

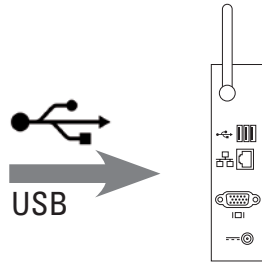

3. Data Transfer Hub

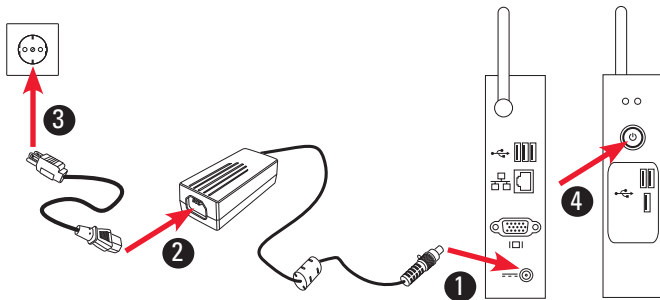

4.



60 sec.

5. App Store



6. WiFi



7. WiFi



8. Leica DMshare



9. Photo

