Nikon Mikroskop / stéréo microscope SM-5

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Prism housings

Looking into the eyepieces, turn the prism housings with both hands symmetrically to bring the right and left circular viewfields into coincidence. (Interpupillary distance adjusting range : 56~75mm)

Eyepieces

10× (Field number 22.7)
Not interchangeable with regular microscope eyepieces.

Diopter ring

Rotate the diopter ring on the left-hand eyepiece, after focusing the specimen through the right-hand by fine focus knobs.

Microscope head lock screw

Release the lock screw to rotate the microscope head horizontally (360° rotatable) or remove it from the focusing mount.

Fine focus knobs

Fine focus range is about 40mm.
For focusing, turn the both knobs in the same direction.
To make the rotation tight, turn the knobs in the direction of the arrow, "TENSION", marked on the knobs. Turning in the opposite direction of the arrow mark slackens the rotation tightness.

Objective

2×, fixed type

Stage plate

Removable by releasing the lock screw. Use its opal or black side according to the specimen to be observed.

Pillar clamp

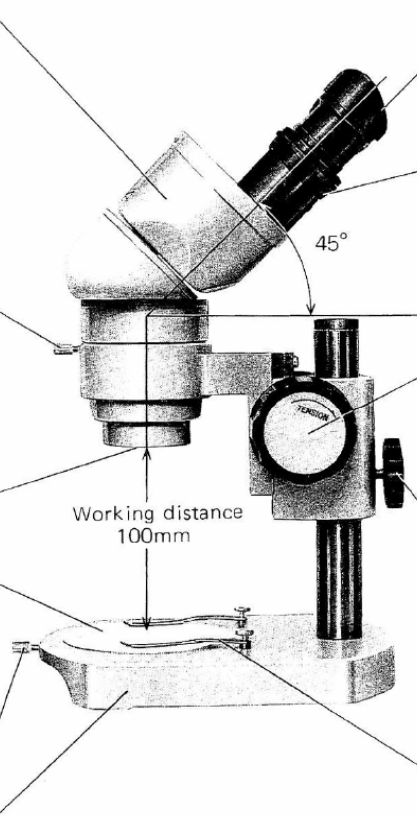
Holding the microscope body with one hand, release the pillar clamp to move the focusing mount along the pillar.
Clamp it tightly in the desired position.

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Ryf AG Grenchen/SO